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## About this document

This document resumes the safety information given in the operating instructions manual for the **Hei-VOLUME Distimatic Pro** automatic module and the translation into all official languages of the European Union.

This document in no way replaces the operating instructions for the product!

For more detailed information, please refer to the original instructions (German language) or the translation of the original instructions (English language) delivered with the product.

## Typographic conventions

Standardized symbols, highlighting elements, and signal words are used in this document to identify warnings, cautions, important information, and special text contents.

Symbol	Signal word / explanatory note
	<p>Warning symbols in combination with a signal word indicate dangers:</p> <p><b>DANGER</b> Indicates a hazardous situation which, if not avoided, will result in death or serious injury.</p> <p><b>WARNING</b> Indicates a possible hazardous situation which, if not avoided, could result in death or serious injury.</p> <p><b>CAUTION</b> Indicates a hazardous situation which, if not avoided, could result in minor or moderate injury or material/environmental damage.</p>
	<p>Mandatory signs are used to indicate important information regarding the product handling.</p> <p>This information is used to ensure operational safety and to maintain the value of the product.</p>

## Copyright protection

This publication is protected by copyright and intended for internal use by the purchaser of the product only.

No part of this publication may be transmitted or reproduced in any form, by any means, without the prior written consent of the copyright owner Heidolph Instruments GmbH & Co. KG. Any violation is subject to compensation for damages.

## Intended use

The **Hei-VOLUME Distimatic Pro** automatic module, according to the EU Machinery Directive 2006/42/EC is a partly completed machine that has been specially developed for the automatic filling and draining of Heidolph rotary evaporators.

The following processes can be controlled with the automatic module:

- filling educt
- drainage of distillate
- drainage of residue (optional)

## General safety information

- Before commissioning and using the device, familiarize yourself with all the safety regulations and guidelines for occupational safety applicable at the place of use and observe them at all times.
- Only operate the device if it is in perfect technical condition. In particular, ensure that there is no visible damage on the device itself and, where necessary, on connected devices or the supply connections.
- If there is missing or misleading information on the device or on occupational safety, contact the responsible safety specialist or our technical service.
- Only use the device in accordance with the regulations for intended use.

## Electrical safety

- Ensure that the voltage indicated on the rating plate matches the supply voltage of the country in which the device is being used.
- Ensure that the mains socket-outlet is protected by means of a residual-current device (RCD).
- Always use the supplied power supply 3-pin cord provided with the device (phase, neutral, ground).
- Prior to use, check that the device and the power supply cord are free of visible damage.
- Have repairs and/or maintenance work on the device carried out exclusively by an authorized and skilled electrician or by the technical service department of Heidolph Instruments.
- Always switch the device OFF and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

## Operational safety

- Do not make any unauthorized changes or modifications to the device!
- Only use genuine spare parts and accessories, or those expressly approved by the manufacturer!
- Rectify malfunctions or faults on the device immediately.
- Shut down the device and disconnect it from the power supply if it is not possible to eliminate the malfunction or rectify the fault immediately.
- Observe all other applicable regulations such as laboratory and workplace guidelines, recognized safety technology rules and special local regulations.
- Observe all relevant general and safety instructions for the connected peripheral devices (observe the supplied documentation!).

## Biohazard

When processing biohazardous substances, take appropriate measures to prevent hazards to persons and the environment, including:

- Instruction of the personnel regarding the necessary safety measures.
- Provision of personal protective equipment (PPE) and instruction of the personnel in its use.
- Marking of the device with the biohazard warning symbol:

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the responsible personnel is the sole responsibility of the operator!

## Work safety

### Assembly

---

#### **CAUTION: Material damage, loss of production**

Due to improper system tubing and wiring as well as unauthorized changes to the system, there is a risk of direct and indirect damage to property!

- For system tubing and wiring, follow all notes and instructions provided in the assembly instructions provided (part number 01-001-009-12-0).
  - Do not make any unauthorized changes to the device, system, tubing, nor the existing cable connections.
  - If necessary, contact our technical service (contact details see operating instructions).
- 

### Control panel

---

#### **CAUTION: Damage to the control panel**

Tip or sharp objects can damage the surface of the display on the control panel!

- To operate the device, touch the display only with your fingertips or a special stylus with a soft-touch tip.
-

## Switch the device on/off

### CAUTION: Property damage, loss of production

If the control box is switched on and off incorrectly, there is a risk that peripheral devices will not be detected correctly or will switch in an uncontrolled manner!

- Observe the order given when switching the control box and the connected peripheral devices on and off.
- In particular, before turning off the control box, make sure that all processes have been completed properly.

Observe the following sequence when turning the automatic module and attached peripherals on and off:

- **TURNING ON THE SYSTEM: First, turn on all peripheral devices** so that they are recognized when initializing the automatic module software.
- **TURNING OFF THE SYSTEM: First, turn off all peripherals** to put the entire system in a safe operating condition.

## Calibrate the flask level sensor

### CAUTION: Loss of production

The function of the level sensor in the evaporation flask (S4) may be affected by electromagnetic fields. This can lead to a loss of the operational quality of the entire system.

- During calibration, commissioning and operation, ensure that there are no sources of interference in the immediate vicinity of the evaporation flask to ensure proper operation of the level sensor at all times.

## System check

### CAUTION: Property damage, loss of production

Faulty connections or signaling can lead to malfunctions or uncontrolled switching operations!

Leaks reduce the performance of the entire system.

- Correct any indicated errors and/or detected malfunctions immediately.
- In the event of a continued fault indication/malfunction, contact your local sales representative or our technical service department (contact details see operating instructions).

## Perform overpressure test

---

### **WARNING: Risk Of Injury, Damage To Property**

Due to faulty components, faulty assembly of components or faulty connections, there is a risk that pressure-applied components will suddenly loosen or rupture during an overpressure test!

- Before performing an overpressure test, ensure that all components are properly installed.
  - In particular, make sure that all screw connections are properly tightened.
  - Before starting the overpressure test, close all safety covers, hoods and doors of the entire system.
  - Ensure that third parties are clear of the danger zone.
  - Use the required personal protective equipment (eye protection, hand protection, etc.).
- 

## Device service

---

### **WARNING: Danger of electric shock**

Live components are installed inside the device.

When opening the device, there is a risk of touching live components.

- Switch the device's main switch off and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

Penetrating liquid poses the danger of an electric shock.

- When cleaning, avoid the penetration of liquids.
- 

## General cleaning instructions

---

### **CAUTION: Damage to the device**

Improper cleaning can damage the surfaces of the device.

Penetrating liquid can damage the electronic components inside the device.

- Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.
  - Never use any aggressive or abrasive cleaning agents or aids.
-

### **Repairs**

Repairs to the device may only be carried out by authorized skilled experts!

Unauthorized repairs during the warranty period will result in the loss of the warranty claim.

The owner is solely liable for damage caused by unauthorized repairs.

In case of repair contact an authorized dealer or our technical service.

### **Maintenance**

The device contains no user-serviceable components. If necessary, in the event of abnormal operating behavior such as excessive noise or heat generation, for example, contact our technical service.

### **Dismantling**

Observe the instructions contained in the assembly instructions supplied when dismantling.

### **Other regulations**

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

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© Heidolph Instruments GmbH & Co. KG